




Instruction Manual



Light Power Meter

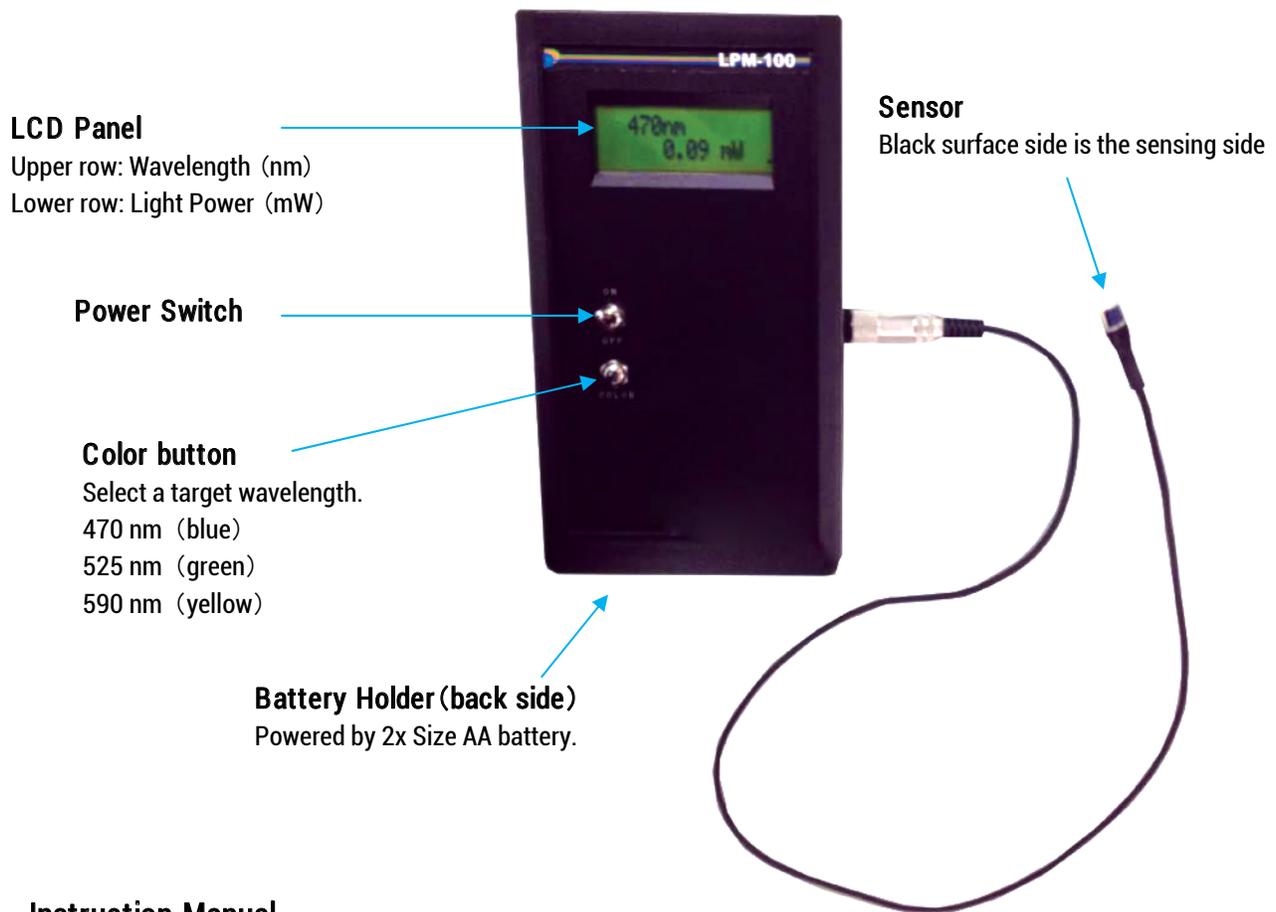

Instruction Manual

- ① Turn on the power switch.
 - ② Push color button several times for setting the target wavelength.
470 nm → 525 nm → 590 nm → 470 nm...
 - ③ Expose the sensor to the target light source. If you measure a light from optic fiber, the fiber end should touch the surface of the sensor at center, vertically.
 - ④ You should see the value of light power from the source on the LCD panel.
- ※ For reducing background of the environmental light, we recommend that the measuring should be done in the box-like cover for making a shade on the sensor.